

KERN & Sohn GmbH

Ziegelei 1 D-72336 Balingen E-Mail: info@kern-sohn.com Phone: +49-[0]7433- 9933-0 Fax: +49-[0]7433-9933-149 Internet: www.kern-sohn.com

Assembly instructions Tripod

KERN PCD-A03

Version 1.0 08/2011 GB





KERN PCD-A03

Version 1.0 08/2011

Assembly instructions Tripod

Scope of delivery:



Qty.	Components
1	Tripod
4	Screws
1	Cable guidance

Technical data

Tripod height	25 cm
Dimensions	195 x 195 x 250 mm

2 PCD-A03-IA-e-1110

Assembly:

There are several possibilities to use the tripod PCD-A03:


Variant 1

The tripod is placed on a base foot next to the balance.

Variant 2:


The tripod is assembled to the platform of the balance

⇒ Slacken 4 screws on the rear side of the display unit and remove the plastic covers [1]

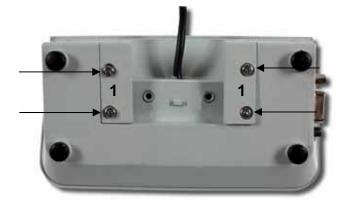


⇒ Plug together display unit and tripod acc. to illustration

Make sure that the cable is passed through the designated opening [2].

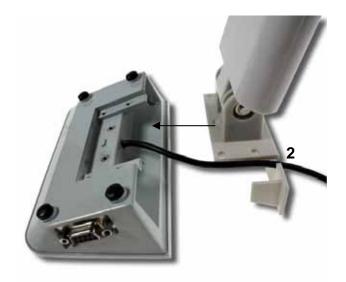


⇒ Tighten 4 screws

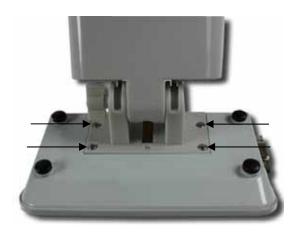

PCD-A03-IA-e-1110 3

Variant 1

⇒ Put the cable on the rear side of the tripod through the cable guidance and fasten cable guidance by 4 screws

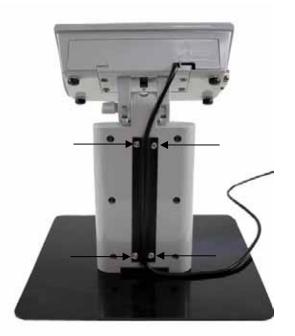



4 PCD-A03-IA-e-1110

Variant 2:

□ Unscrew 2 screws at the lower side of the adjustment foot



⇒ Push together tripod and support of the platform



PCD-A03-IA-e-1110 5

⇒ Fasten with screws



⇒ Put the cable on the rear side of the tripod through the cable guidance and fasten cable guidance by 4 screws



6 PCD-A03-IA-e-1110



PCD-A03-IA-e-1110 7